Levenhuk 2S/3S biological microscope

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levenhuk[®]
Zoom&Joy

- 1 Monocular head
- 2 Microscope tube
- 3 Objective lens
- 4 Stage
- 5 Disc diaphragm
- 6 Mirror
- 7 Eyepiece
- 8 Focus knob
- 9 Microscope arm
- 10 Stage clips
- 11 Microscope base

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Congratulations on your purchase of this wonderful Levenhuk microscope!

This instrument was made to the highest international standard and, when used in accordance with this manual, is not hazardous to the environment, your possessions or your health.

This microscope was designed for observations of biological specimens and for basic lab work.

Specifications

Magnification	200x	Focusing range	40 mm (1.58 in)
Achromatic objective lens	6x	Stage size	90 mm x 90 mm (3.54 x 3.54 in)
Eyepiece	2.5x	Mirror diameter	40 mm (1.58 in)
Working distance	4.5 mm (0.18 in)		

Using the microscope

- Place your microscope on an even surface. The mirror acts as an illumination source, so make sure that you choose your working place accordingly. During daytime an ideal spot is by the window, in the evenings it should be well-lit by a bright lamp. Rotate the mirror, so that you see a bright circle of light through the eyepiece. Do not, under any circumstances, point the mirror at the sun, as this may cause permanent retinal damage.
- Place a prepared slide on the stage and secure it in place with stage clips.
- Use the focus knob until the objective lens almost touches the cover glass of the slide.
- Avoid touching the slide with the objective lens, as this may damage on or the other. After lowering the microscope tube, look through the eyepiece, then use the focus knob to bring the tube up and focus the view.
- For additional comfort during your observations you can incline the eyepiece and the stage by pulling the microscope arm.

Care and maintenance

- After unpacking your microscope and before using it for the first time check for integrity and durability of every component and connection.
- Do not touch the optical surfaces with your fingers. You can use a soft napkin or a cleaning wipe, dipped in ether or absolute alcohol, to clean lenses. We recommend using special optics cleaning fluids from Levenhuk. Do not use any corrosive or acetone-based fluids to clean the optics.
- Abrasive particles, such as sand, should not be wiped off lenses, but instead blown off or brushed away with a soft brush.
- Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws.
- Protect the device from sudden impacts and excessive mechanical force.
- Store the device in a dry, cool place away from dust, hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures. Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity.
- Be careful during your observations, always replace the dust cover after you are finished with observations to protect it from dust and stains.
- If you are not using your microscope for extended periods of time, store the objective lenses and eyepieces separately from the microscope.

International Warranty

- The Vendor guarantees that the quality of the Levenhuk product purchased complies with the technical documentation requirements on conditions that the consumer observes rules of transportation, storage conditions and operating instructions.
- Levenhuk Ltd. warrants the product against defects in materials.
- During the warranty period the Buyer can return the defect product to the Vendor or to Levenhuk service center. Levenhuk Ltd. will repair or replace the defect product at its own discretion.
- No claims are accepted in case the properly filled warranty slip is absent or contains corrections, or if the defect product has not been provided by the Buyer. The manufacturer or the seller is not liable for any damages caused by misuse of the product. No guarantee is provided in cases the product was used for purposes other than that intended or has mechanical damages, scratches, cracks, optics damages, as well as in cases the product is broken-down due to impact damages, squeezing or stretching or the product has been disassembled or repaired by unauthorized personnel.
- This warranty does not cover components and/or accessories with limited use period, batteries, etc.

The period of warranty is 3 (three) years beginning on the date of purchase. Please keep the warranty slip along with your receipt.

For any warranty services please contact the local Levenhuk branch.

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